
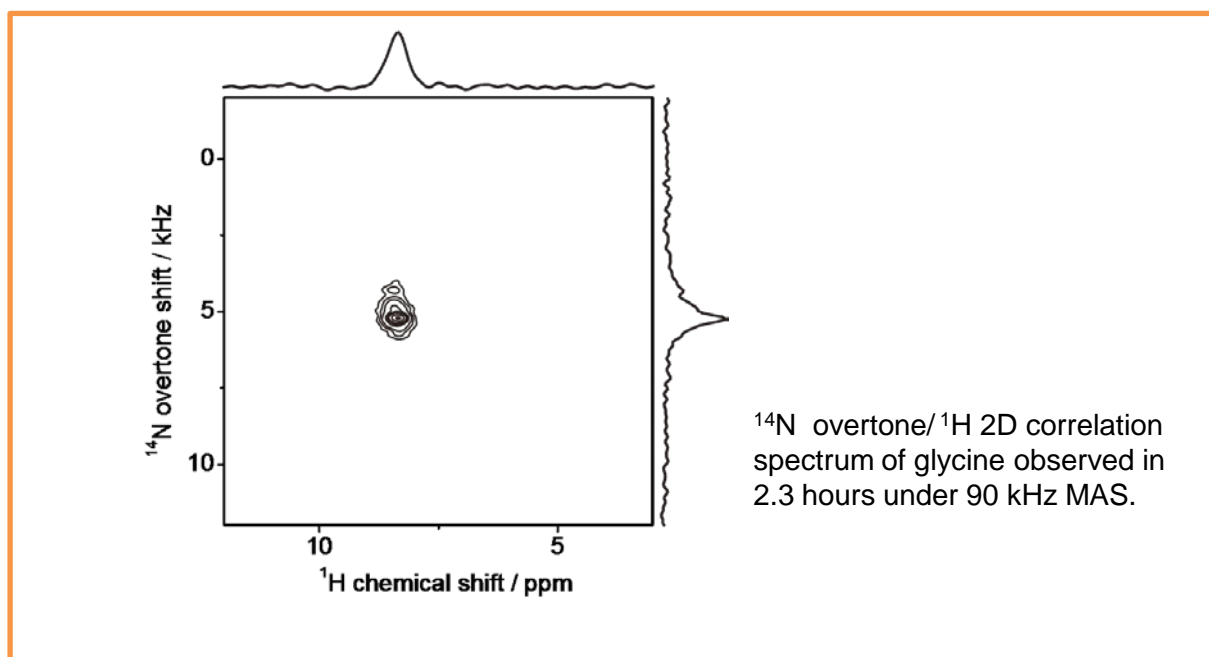


Overtone solid-state NMR spectroscopy on Nitroge-14

Product used : Nuclear Magnetic Resonance Spectrometer (NMR)

^{15}N NMR is widely used because of the importance of nitrogen in chemistry, materials, biology, environment etc. However, very low abundance of ^{15}N (<0.4%) results in poor sensitivity and thus makes observation time-consuming. On the other hand, the rest of nitrogen atoms are also NMR sensitive nucleus of ^{14}N . Despite the high abundance of ^{14}N (>99%), it's application is rather limited due to the huge quadrupolar interactions and its spin quantum number $I = 1$. The introduction of very fast MAS (>70 kHz) enabled the ^1H detected $^{14}\text{N}/^1\text{H}$ correlation spectroscopy (Nishiyama et al., JMR 208 (2011) 44-48). However, it demands precise magic-angle adjustment, moreover, the sensitivity is largely affected by the molecular motion. It was shown that the use of ^{14}N overtone transition avoids these difficulties under MAS (O'Dell et al. Chem. Phys. Lett. 514 (2011) 168-173) with a cost of sensitivities. Since the overtone transition between -1 and +1 energy level is forbidden, the low sensitivity fundamentally comes from small transition moment between these levels. Here, we have developed ^1H detected ^{14}N overtone/ ^1H correlation spectroscopy under ultrafast MAS conditions > 70 kHz to achieve highly sensitive ^{14}N NMR spectroscopy of rigid solid.

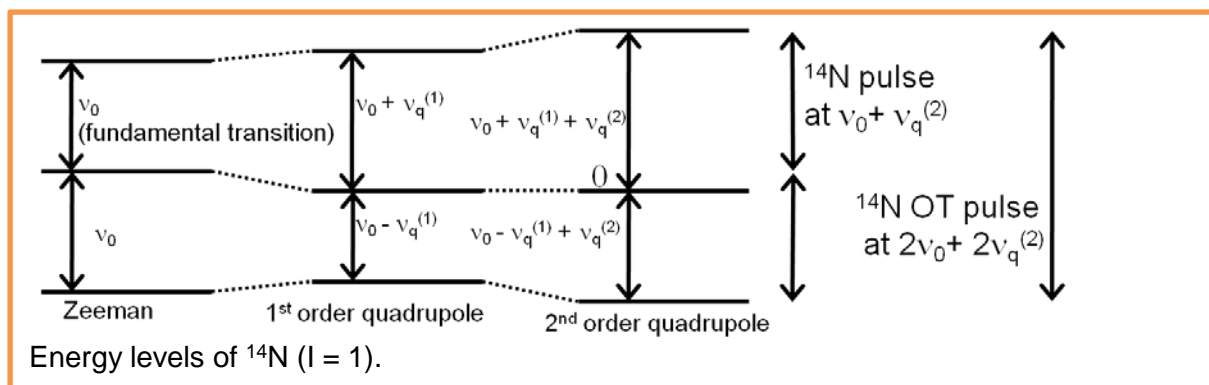

Y. Nishiyama, M. Malon, Z. Gan, Y. Endo, T. Nemoto, J. Magn. Reson. 230 (2013) 160-164.

Copyright © 2018 JEOL Ltd.

Certain products in this brochure are controlled under the "Foreign Exchange and Foreign Trade Law" of Japan in compliance with international security export control. JEOL Ltd. must provide the Japanese Government with "End-user's Statement of Assurance" and "End-use Certificate" in order to obtain the export license needed for export from Japan. If the product to be exported is in this category, the end user will be asked to fill in these certificate forms.

